

# Palintest

Water Analysis Technologies

## Ozone

Photometric test for Ozone in drinking and treated water

**COLOUR CHANGE**

Colourless to Pink

**RANGE:**

0 - 3 mg/L O<sub>3</sub>

**TECHNICAL RESOURCE:**

Ozone Test Method, Technical Information

01



Fill to the 10 mL line with sample.

02



Use this to "blank" the instrument. Place in cell holder and press Blank.

03



Pour out sample leaving a few drops.

04



Add a DPD 4 tablet to those drops. Then, crush to form a paste.

05



Fill to the 10 mL line with further sample and stir.

06



Place the tube in the cell holder and press Measure to obtain the Ozone concentration.